

Instructions for use


Rotor F-45-72-8, Rotor F-45-48-11

English (EN)

This document complements the operating manual for the Concentrator/Vacufuge® and does not replace it. Therefore, please also read the operating manual for the Concentrator before starting up the rotors for the first time.

1 Using the rotor feet

The rotor feet prevent the vessels from sliding out of the bores when positioning the rotor. The rotor feet may remain in place during centrifugation.

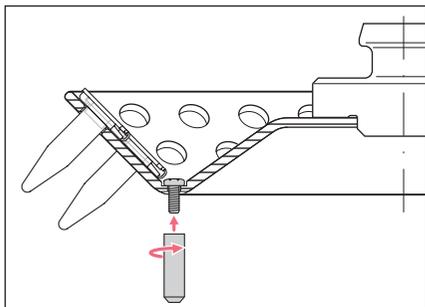

- ▶ Insert the 3 screws through the bores.
- ▶ Tighten the rotor feet.

2 Using the spacer

Use the spacer to centrifuge with 2 rotors at the same time. The rotors can be positioned freely. The following maximum loads apply:

- Lower rotor: $\frac{2}{3}$ of the max. load
- Upper rotor: $\frac{1}{3}$ of the max. load

3 Ordering information

Order no. (International)	Description
5490 034.007	Rotor F-45-72-8 72 positions for 0.5 mL tubes
5490 030.001	Rotor F-45-48-11 48 positions for 1.5/2.0 mL tubes
5301 316.005	Spacer for simultaneous operation of two rotors for F-45-72-8 and F-45-48-11
5490 030.800	Rotor feet for F-45-72-8 and F-45-48-11 Set of 3